

Introducing a truly integrated Flow Cytometry Reporting solution for Anatomic Pathology laboratories

The days of transcribing individual marker results into the pathology report and passing around histograms and scatter plots are over! NovoPath™, a leading AP Software Platform, now offers a simple and direct way to retrieve and parse the results by defining specific panels with any combination of markers that mirror the results being downloaded from the instrument.

- Provides complete setup of individual markers and their inclusion and grouping into standard run panels for specific disease entities for maximum efficiency.
- Allows users to add new gates per panel as needed and to order add-on panels to provide additional information to the diagnostician.
- Provides automatic download of all histogram and dot plot images that will be attached directly to the patient's case so that they are retrievable for viewing, printing and inclusion in the patient report if desired.
- Predefined dictionary results can be setup and made available for quick diagnosis entry.
- Calculate QC Result data based on the marker data received from the instrument to save technologist time.
- Custom define patient report templates to display flow data as desired for a better competitive advantage in the marketplace.

Analysis [Case #: FC09-00149/Status: Provisionally Diagnosed]

Patient: Peter M. Anderson

Client: Viability Gated C

HAIRY CELL Panel

Regate Gate # 1 Lab Completed

PDF Add to Report

Amendment/Addendum --SELECT--

Marker	Value	Status
CD2	12	L
CD3	45	L
CD4	11	L
CD5	5	
CD7	65	
CD8	20	
CD10	44	
CD13	12	
CD14	31	
CD19	78	H
CD20	45	M
CD23	45	
CD33	45	
CD34	45	H
Cd38	32	
Cd45	34	
CD56	65	
HLA-DR	3	
FMC7	89	
Kappa		
Lamda		
CD20/Cd38	37	
CD25		

Analyzed Region	Result	Range
Analyzed Region		NONE
CD4+CD8=CD3	68.889	85-115
CD2=CD3	0.267	90-110
T+B+NK+BLASTS	188.000	85-115
CD19=CD20	173.333	85-115
CD45 from Gate B	34.000	50-100

CONTACT:

To find out more about the Flow Cytometry Module, contact your Sales Representative at: 877-668-6123 or email us at: sales@novopath.com

You can include all flow, cytogenetics and biopsy results on the same report, offering true multi-modality reporting!